**Informationsbroschüre des Herstellers  
Information supplied by the manufacturer**

## **EKASTU-Augendusche Aqua Push EKASTU-Eye Wash Unit Aqua Push (Best.-Nr. / Order no.: 177 500)**

### **DIN EN 15154-2**

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**The EKASTU-Eye Wash Unit Aqua Push complies with DIN EN 15154-2 and the BGI/GUV-I 850-0 "Safe working in laboratories - principles and guidelines" issued by the German Social Accident Insurance (DGUV). Please observe the requirements of this standard and guideline as well as any other applicable regulations.**

**EU-Declaration of Conformity is available for download from [www.ekastu.de](http://www.ekastu.de).**

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## 1. General notes

The use of the EKASTU Aqua Push eye shower requires full knowledge of and compliance with this information brochure. The EKASTU Aqua Push eye wash is intended exclusively for the purposes stated in the information brochure. These is used to rinse out foreign substances from the eye as part of first aid treatment. It does not replace an ophthalmological examination, but can help to minimize injuries. We therefore recommend seeking medical help after an eye accident.

EKASTU Safety GmbH assumes liability within the scope of its general terms and conditions. If you do not have them, they will be sent to you on request. The manufacturer is not liable for damage caused by non-compliance with the information brochure.

## 2. Use, Storage, Maintenance

Eye showers with a water connection are mandatory first aid equipment for workplaces where hazardous substances are handled. The eye shower is used for targeted rinsing of the eyes and for rinsing the face.

Store the EKASTU Aqua Push eye shower and spare parts in a dry place, protected from dust and corrosive and oxidant agents.

The EKASTU Aqua Push eye shower must be maintained in accordance with the following maintenance schedule:

Type of inspection	Time limits	Implementation by User/Instructor/Specialist	Type of work to be carried out
Function test of the device	weekly	User/ Instructor/ Specialist	<ul style="list-style-type: none"><li>• Check for contamination and leaks</li><li>• Check shower head for calcification</li><li>• Check the water supply hose for twists or kinks</li><li>• Operate the eye shower for approx. 5-10 seconds <sup>1)</sup></li></ul>
Maintenance of complete equipment	yearly	Instructor/ Specialist	<ul style="list-style-type: none"><li>• Check for dirt (e.g. dust or calcification deposits), leaks, closing behavior of the release handle</li><li>• Check the water supply hose for twists, kinks or other damaged areas</li><li>• Replace defective parts</li><li>• Operate the eye shower for approx. 2 minutes <sup>2)</sup></li></ul>

<sup>1)</sup> This will check the function and replace the stagnant water in the eye shower with fresh water and thus prevent germination.

<sup>2)</sup> This prevents the standing water from germinating.

## 3. Installation site and Marking

The installation location of the eye shower should be chosen carefully, because every second counts in the event of an eye accident, so it is crucial that the shower is at least in the same room and that the distance from the workplace is no more than 10 meters or less or that the eye shower can be reached within 10 seconds or less. The location of the eye shower must be known to all employees. Access must be kept clear at all times and a free space of at least 15 cm must be maintained around the eye shower. In addition, the installation location of the eye shower should be marked with a safety sign. To do this, use the enclosed self-adhesive sign 150 x 150 mm and attach it at eye level. In order to avoid damage or consequential damage caused by leaking water, the eye shower should ideally be installed in the area of a sink; if this is not possible, the water can be safely collected in drip trays during the rinsing process.

## 4. Connection and Installation

The eye shower should be connected to a cold drinking water connection. Connecting the eye shower to a cold drinking water connection prevents burning with excessively hot water and/or the formation of germs, and the cold water also has a pain-relieving effect in the event of burning and/or burns to the eyes.

To ensure proper function, the flow pressure must be at least 4.5 bar and no more than 5.5 bar.

## 5. Mounting

Wall mounting is carried out on a suitable surface (wall) using a wall bracket supplied.

The required fixing technology must be specified by the installer depending on the construction.

## 6. Malfunction

The eye shower must be kept in a functional condition. In the event of a malfunction of the eye shower, the responsible person must be informed and the necessary repair or replacement of the eye shower must be initiated. Employees must be informed for the duration of the repair or replacement of the eye shower.

## 7. Disposal

The eye shower must be disposed of in accordance with the applicable local waste disposal regulations.

Informationsbroschüre\_ EKASTU-Augendusche Aqua Push\_177 500\_DE\_EN\_01-06.25

**EKASTU Safety GmbH (Group)**  
Schänzle 8 • D-71332 Waiblingen  
Germany

**Tel.: +49(0)7151 975099-0**  
**Fax: +49(0)7151 975099-30**

**Internet: [www.ekastu.de](http://www.ekastu.de)**  
**eMail: [info@ekastu.de](mailto:info@ekastu.de)**

**EKASTU Safety AG**  
Kreuzbühl 8 • FL-9493 Mauren  
Principality of Liechtenstein

**Internet: [www.ekastu.li](http://www.ekastu.li)**  
**eMail: [info@ekastu.li](mailto:info@ekastu.li)**  
**Tel.: +423 380 06 06**